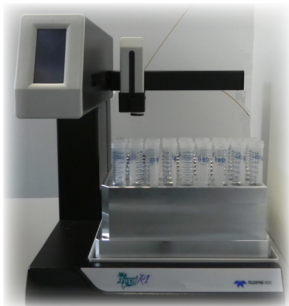
Foxy R1 Cooling

- ▶ Beiblatt zur Installation der Kühloption
Supplement Installing Cooling Option

V7034

HPLC

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Table of Contents

Note **For your own safety, read the manual and always observe the warnings and safety information on the device and in the manual!**

German Table of Contents	4
Intended Use	20
Overview of Foxy R1 Cooling	21
Safety	22
Symbols and Signs	22
Scope of Delivery	23
Checking the Scope of Delivery	23
Installation	24
Installing the Cooling Option	24
Mounting the Cooling Plate	24
Inserting the Rack	27
Assembling/Disassembling the Hood	30
Replacing the Rack	30
Troubleshooting	31
Possible Problems and Rectifications	31
Further Measures	31
Contacting the Technical Support	31
Technical Data	32
Requirements for the Thermostat	32
Accessories and Spare Parts	33
Abbreviations and Terminology	34

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Intended Use

The KNAUER cooling option for fraction collector Foxy R1 serves for the cooling of fraction samples.

You need an external thermostat to transport the coolant at 4 °C into the cooling plate.

The cooling option consists of three subassemblies:

- The cooling plate with connection for the coolant
- Optionally, different racks for the fractions are available:
 - Rack for Eppendorf vials (12 × 12)
 - Rack for Falcon tubes (8 × 9)
 - Cooling adapter for 96 well plates
- The hood as cooling cover to prevent the condensation of water and to cool the fraction sample at a steady temperature

Necessary tools for an installation:

- Scissors to shorten cooling tubes and insulation tubes
- Pliers to secure the cooling tube to the olive-type tube with cable binder
- Coolant (mixture of water and antifreeze) to avoid ice formation inside the device.

Overview of Foxy R1 Cooling

Fig. 7 Example for cooling after removing the hood

Legend

- ① Diverter valve assembly, mechanically adjustable height
- ② Fraction tubes
- ③ Rack for Falcon tubes
- ④ Cooling plate with connectors for cooling tubes

Safety

Note **Keep this supplement on the installation of the cooling option and store it together with the user manual for the fraction collector Foxy R1.**

When installing or replacing the heavy racks or the cooling plate, pay attention to the safety notes.

Professional group

Always have the KNAUER cooling option for the fraction collector installed and operated by trained technicians, e.g. chemical engineering assistants.

If you do not belong to this or a comparable professional group and do not fulfill these requirements, under no circumstances may you perform the work described in this service supplement. In this case, please contact your superior.

Symbols and Signs

There are no warning or other signs to be seen on the cooling of the fraction collector.

Warning signs in the supplement	Explanations
	Hand injuries during installation possible

Scope of Delivery

Note: Only use original parts and accessories made by KNAUER or a company authorized by KNAUER.

Delivery	Hood	<input type="checkbox"/>
	Cooling plate	<input type="checkbox"/>
	Rack	<input type="checkbox"/>
	Service supplement	<input type="checkbox"/>
Accessories kit	Accessories kit for cooling tubes	<input type="checkbox"/>

Checking the Scope of Delivery

1. Check whether the device and accessories are complete.
2. If a part is missing, inform the technical support of KNAUER (see page 31).

Installation

Installing the Cooling Option

Prerequisite

- The device has been switched off.
- The power plug and the LAN cable have been removed.
- The diverter valve assembly has been brought to the upmost position.

Mounting the Cooling Plate

Process

1. Place the cooling plate ④ onto a table.
2. The connectors for the cooling unit point to the back.
3. Using scissors, cut the cooling tube ② to the required length.
4. Using scissors, cut the insulation tube ① to the required length and push it onto the cooling tube.
5. Attach the cooling tube to the olive-type tube ③ of the cooling plate.
6. Accordingly attach the second cooling tube and the insulation tube.
7. Using pliers and cable binders, secure the cooling tubes to the olive-type tubes ③.

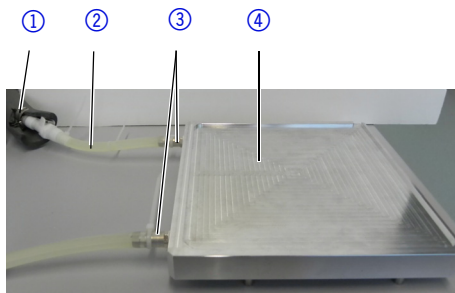
Process

Fig. 8 Cooling plate and rear side connectors for cooling

8.  **Caution!**

Hand injuries are possible.

Consider the high weight of the cooling plate. Place the cooling plate onto the drip tray, consider the orientation of the rack-guide pins ⑤.

Process

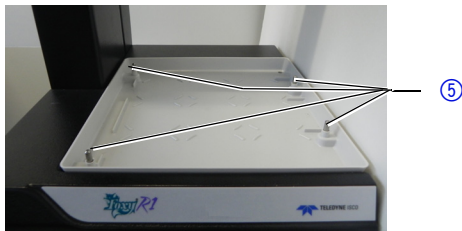
Fig. 9 Foxy R1 drip tray with rack-guide pins

Inserting the Rack

Note **The rectangular corner of the rack must be in the front on the right hand side. Otherwise, the rack cannot be mounted correctly.**

The rack lies flat on top of the cooling plate and is not attached to it.

Process

1.

Caution!

Hand injuries are possible.

Consider the high weight of the rack. Start with positioning three-quarters of the rack on the cooling plate. Then push the rack manually further onto the cooling plate until cooling plate and rack ① fit perfectly.

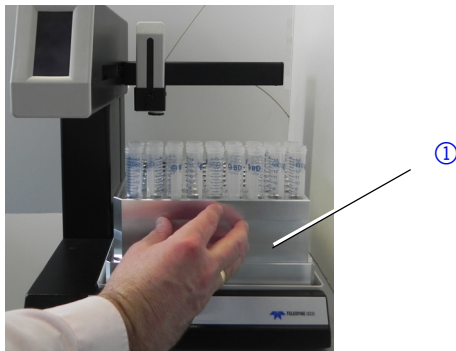
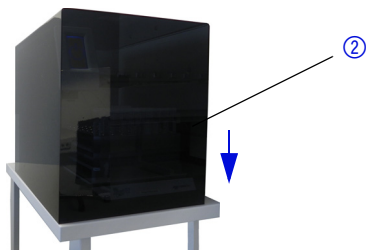
Process

Fig. 10 Pushing the rack into its end position

2. Adjust the height of the diverter valve assembly accordingly to the height of the sample vials.
3. Set on the hood ② vertically from top down, the holes for the cooling tubes point to the back (see arrow). The hood is not screwed to the fraction collector.

Process**Fig. 11** Hood on top of the rack

Assembling/Disassembling the Hood

The hood helps to cool the fraction samples and prevents the development of condensed water.

Always assemble the hood vertically from top down with the holes for the cables and cooling tubes pointing to the back.

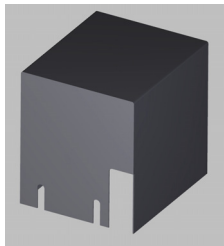
Fig. 12 Hood rear side with holes for cables and cooling tubes
To disassemble the hood, lift it vertically from bottom up (see arrow).

Note Consider that you have enough space for the assembly/disassembly.

Replacing the Rack

Lift the rack with both hands and pull it towards the front. Then grip the rack with both hands and remove it.

Troubleshooting

Possible Problems and Rectifications

Problem	Solution
Ice formation in thermostat	Raise the amount of antifreeze.
Insufficient cooling	Raise the circulation velocity of the thermostat.

Further Measures

1. Check all cabling.
2. Check all screw fittings.
3. Check device for leaks.

Contacting the Technical Support

You have various options to contact the Technical Support:

Telephone +49 30 809727-0

Fax +49 30 8015010

E-Mail support@knauer.net

You can make your requests in English and German.

Technical Data

Weight (with cooling option, without racks)	12.0 kg
Dimensions (width x height x depth)	15 cm x 10 cm x 15 cm
Space requirements (width x height x depth)	35 cm x 50 cm x 50 cm
Temperature range	2°C - 60°C
Materials	Anodized aluminum

Requirements for the Thermostat

Cooling capacity	2 °C
Flow rate (suction)	20 liters/minute

Accessories and Spare Parts

For repeat orders of spare parts use the enclosed packing list. Contact the Technical Support in case there are any questions on spare parts or accessories.

Name	Order number
Hood, complete	A59115
Cooling plate, complete	A59116
Rack with cooling option (12 x 12)	A59117
Rack (8 x 9)	A59118
Cooling adapter Foxy R1	A59119
Supplement DE/EN cooling option	V7034
Accessory kit for Foxy R1	F59100-1
Vinyl tubing for funnel rack	A59120
Diverter valve	A59121

Abbreviations and Terminology

Term	Explanation
Tube/Vial/Well	Containers of different sizes and shapes

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© Wissenschaftliche Gerätebau
Dr. Ing. Herbert Knauer GmbH
All rights reserved.
The information in this document is
subject to change without prior notice.
Translation of the original German edi-
tion of this supplement, version 1.0.
2013-04-16
Printed in Germany on environmentally
friendly paper from sustainable forests.

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